

OBIS™ LX/LS and LX SF Quick Start Guide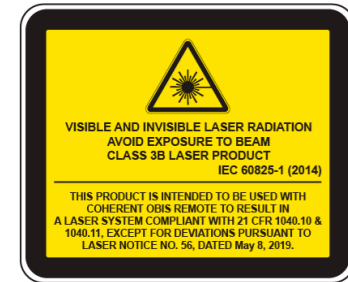

Laser Safety

Laser light, because of its special properties, has safety hazards not associated with light from standard sources. The safe operation of lasers requires that all laser users, and everyone near the laser system, know the dangers involved. The safe use of the laser needs the user to be familiar with the instrument and the properties of coherent, intense beams of light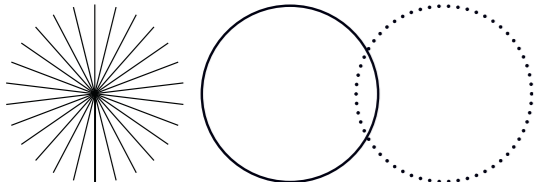

DANGER!
Direct eye contact with the output beam from the laser causes damage and possible blindness.

Laser beams can cause ignition of volatile material such as alcohol, gasoline, ether and other solvents, and can cause damage to light-sensitive material in video cameras, photomultipliers and photodiodes.

Reflected beams can also cause damage.

Laser Safety Precautions

These **safety precautions** are recommended:

- Obey all safety precautions in the operator's manual.
- Caution must be used when solvents are used in the area of the laser.
- Limit access to the laser to approved users who know laser safety procedures and who do not know the risk involved.
- Do not look directly into the laser light source or at scattered laser light from a reflective surface. Do not look down the beam into the source.
- Maintain experimental setups at low heights to prevent accidental beam to eye contact at eye level.
- As a precaution against accidental exposure to the output beam or its reflection, individuals that operate the system must wear laser safety glasses.

WARNING!
While laser safety glasses give protection to eyes from potential damage from exposure, glasses may block light at the laser wavelengths, and that prevents the operator from being able to see the beam. Use strong caution even while safety glasses are worn.

- Use the laser in a closed room. Laser light stays collimated over long distances and shows a possible hazard if not confined.
- Post warning signs in the area of the laser beam to give caution to individuals in the area.
- Recommend to all individuals that operate the laser of these precautions. It is good procedure to operate the laser in a room with controlled and restricted access.

Electrical Safety

The OBIS laser does not contain hazardous voltages.

Do not disassemble the enclosure. There are no user-serviceable components. All units are made to be operated as assembled. The warranty will be voided if the enclosure is disassembled.

Electrostatic charges as high as 4000 volts easily collect on the human body and equipment when handling a device during installation or use and can discharge without detection.

CAUTION!
Take correct ESD precautions during installation and operation to prevent damage or performance degradation.

The laser has input protection. However, permanent damage can occur on devices that receive high-energy electrostatic discharges.

OBIS Laser Installation

This section shows how to set up the OBIS laser to run in Continuous Wave (CW) mode.

The user is responsible for compliance with all applicable laser safety requirements.

For more information about modulation, interface, installation, heatsink use, packaging, and accessories, see the *OBIS LX/LS Operator's Manual* (P/N 2283534).

The OBIS automatically starts laser emission after DC power is supplied. The user is responsible for compliance with all applicable laser safety requirements.

Install Coherent Connection Software

Install the software on a Windows (64-bit) host PC or laptop before the laser is set up and powered on.

- Download the *Coherent Connection* software (which also contains all the necessary drivers) from the Coherent website:
<https://www.coherent.com/resources>

- Close all programs.
- On the PC, double-click this file to start the installation process:
`Coherent_Connection_Setup_v6.0.x.xx.exe`
- Follow the on-screen instructions to complete software installation.

For more software installation and operation instructions, refer to the *Coherent OBIS LX/LS Operator's Manual* (P/N 2283534), available in PDF format on the Coherent website: <https://www.coherent.com/resources>

Required Materials

An optical plate with mount holes at a ~1" distance apart **is not** a correct mount surface without an OBIS Heat Sink. See the *OBIS LX/LS Operator's Manual* (P/N 2283534) for full instructions on heat sink requirements and procedures to install the laser.

A USB cable is also necessary to connect the OBIS laser to a host PC or laptop.


Attach Laser to Heatsink

WARNING!
DO NOT use thermal grease or thermal compounds. The use of thermal grease or compounds voids the Coherent warranty.

- Use a 2.5 mm hex wrench and tighten the screws in a diagonal pattern onto the heat sink.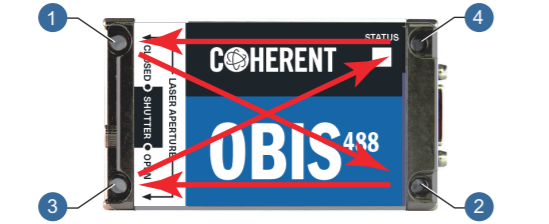

- Torque the mounting screws, initially, to 0.25 N·m (35.4 oz·in.) in a 1-2-3-4 sequence. Use the same diagonal pattern for the last torque setting of 1 N·m (141.6 oz·in.).

Option 1: Connect Laser Without Remote


Attach to Power Supply Direct for use **without** remote:

- Remove the yellow 'OEM Use Only' label from the back of the OBIS laser to access 12V OEM power supply connector. See figure.


CAUTION!
The OBIS automatically starts laser emission after DC power is supplied.

- Attach the USB cable from the host computer to the USB connector on the laser for use **without** remote.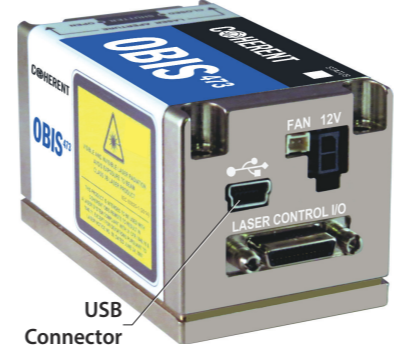

- Insert the 2-pin 12 VDC OEM Power Supply connector into the power connector on the back panel of the laser.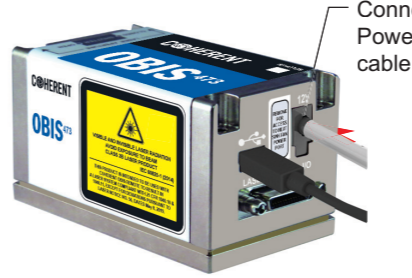

Option 2: Connect Laser With Remote

Perform remote back panel connections.


- Connect the power supply to the remote.

- Attach the SDR cable to the remote.

- Attach other end of SDR connector directly into SDR connector on the laser for use **with** remote: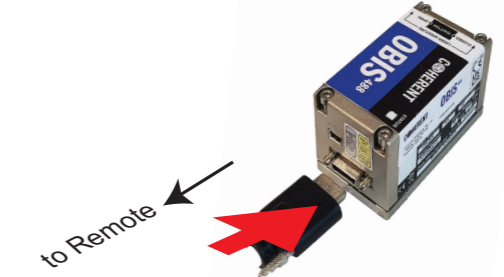

- Tighten all connector thumbscrews. (continue to page 14)

部件名称 Part Name	有害物质 Hazardous Substances					20 Recycling symbol
	铅 (Pb)	汞 (Hg)	镉 (Cd)	六价铬 (Cr(VI))	多溴联苯 (PBB)	
印刷电路板组装 PCB Assembly	X	O	O	O	O	O

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
Perform remote front panel connections:

- Put the interlock connector into the remote.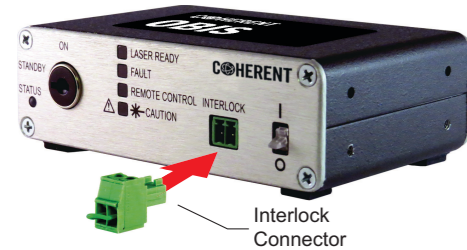

- Make sure that the key is inserted and set the keyswitch to the ON position.

Laser Operation

OBIS: Slide the lever right (to OPEN position) to open the mechanical front shutter to allow light output.

OBIS Fiber Pigtail (FP): Remove the fiber shutter cap.

Shutter Cap removed

To automatically start emission:

- Give 12 VDC power to the OBIS.

Auto Start is enabled as a factory default setting and laser emission happens when power is applied to the laser and thermal stability is achieved. Note that when Auto Start is enabled, there is no 5-second delay.

OBIS starts to emit light after it is thermally stabilized. During stabilization, the LED indicator on top of the laser flashes green.

LED Indicator

When stabilized, the LED indicator changes to white and laser emission starts. Light emission starts at the last power set point.

If emission does not start within 5 minutes and the LED indicator on top displays in solid blue, the Auto Start feature is likely not in enabled. This means the laser has stabilized thermally but is waiting for a command to start emission.


To manually start emission if emission does not autostart:

- 1) Connect the OBIS laser to a PC. Use a USB Type A-to-USB Type Mini B cable.
- 2) Make sure that the OBIS is correctly connected to its 12 VDC power supply.
- 3) Start the Coherent Connection software.
- 4) Press the **Start** button in the Main tab.
- 5) To enable the Laser Auto Start feature, go to the Advanced tab in Coherent Connection and put a check mark in the **Auto Start** box.

Configure Wavelength Settings (OBIS LX Single Frequency Lasers Only)

Use the Operating Power tab in the software to fine-tune the laser's wavelength:

- Adjust the diode temperature.
 - Adjust the operating power to adjust the diode current.
- or
- A combination of both.

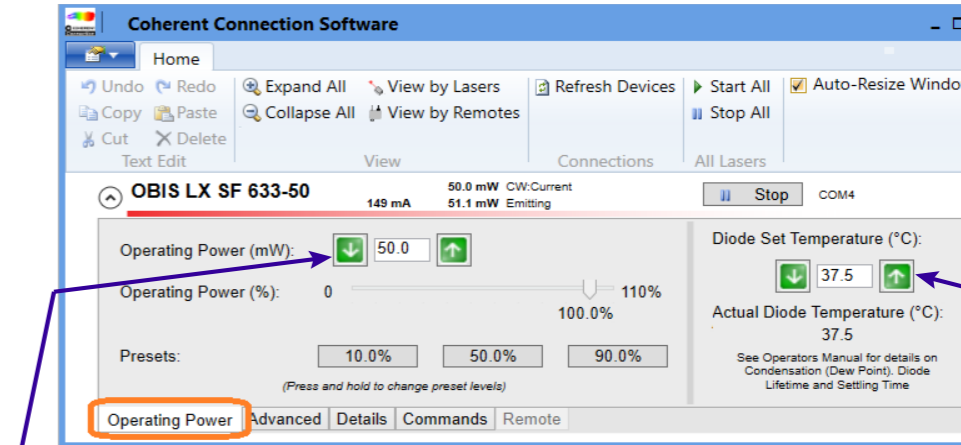Arrows change power level by 1mW (or) Enter power level manually

Arrows change diode temp by 0.1°C (or) Enter diode temperature manually from 20°C to 40°C

The screen below shows the status of the laser:

NOTICE

For Single Frequency Laser, make sure that the laser operates in 'CW:Current' (constant current) mode.

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© Coherent Corp., September 2023 (RoHS)